

Enriched Seawater Medium Recipe

Directions

Modification of L. Provasoli's ES-enrichment for seawater a.k.a. PES (Bold & Wynne 1978). General purpose marine medium for xenic cultures.

1. Aseptically add 20 mL of sterile ES Enrichment Solution per liter of Pasteurized Seawater (30 ppt).
2. Store at refrigerator temperature.

#	Component	Amount	Stock Solution Concentration	Final Concentration
1	Pasteurized Seawater	1 L		
2	Enrichment Solution for Seawater Medium	20 mL/L		