

ES/2 Enriched Seawater Recipe

Directions

Suitable for some of J.A. West's tropical green isolates that require less of the ES-enrichment in order to maintain the desired morphology and rate of growth (West 1984, pers. Comm.).

1. Aseptically add 10 mL of sterile ES Enrichment Solution per liter of Pasteurized Seawater (30 ppt).
2. Store at refrigerator temperature.

#	Component	Amount	Stock Solution Concentration	Final Concentration
1	Pasteurized Seawater	1 L		
2	Enrichment Solution for Seawater Medium	10 mL/L		