

Liver extract infusion Recipe

Directions

For 50 mL Total

1. To approximately 40 mL of dH₂O, add 2.5 g of BACTO LIVER while stirring continuously.
2. Bring the total volume to 50 mL.
3. Cover the flask, and place it into a 50° C water bath for approximately 1 hour.
4. Transfer the flask into an 80° C centigrade water bath for 3-5 minutes.
5. Filter extract.

#	Component	Amount	Stock Solution Concentration	Final Concentration
1	Liver extract (Difco Bacto 0133)	2.5 g/50 mL		