

Malt Medium Recipe

Directions

Suitable for all Prototheca cultures.

For 500 mL Total

1. To 500 mL of dH₂O add the following component.

*For 1.5% agar medium add 7.5 g of agar into the flask; do not mix.

2. Cover and autoclave medium for no longer than 15 minutes to avoid caramelizing sugars.

3. Store at refrigerator temperature.

#	Component	Amount	Stock Solution Concentration	Final Concentration
1	Malt extract (Sigma M-0383)	33.6 g/500 mL dH ₂ O		