

## 2/3 Enriched Seawater Medium

For marine algae that require salinities at **20 ppt**, the basic Enriched Seawater Medium is combined with sterile distilled water prior to inoculating the new culture, such that 2/3 Enriched Seawater Medium combines Enriched Seawater Medium and sterile distilled water **2:1**.

### Directions

#### For 1 L Total

1. Combine components.
2. Store at refrigerator temperature.

#	Component	Amount	Stock Solution Concentration	Final Concentration
1	Enriched Seawater Medium	665 mL		
2	sterile dH <sub>2</sub> O	335 mL		